# SYNTHETIC PYRETHRINS/C-ELISA

## Contents

----

.....

A. Introduction
B. Equipment 2
C. Reagents 3
D. Standards
E. Extraction Procedure
F. Analytical Quantitation9
G. Calculations
H. Hazard Analysis 17
I. [Reserved]
J. Quality Assurance Plan

#### A. INTRODUCTION

1. Theory and Structure

Rendered fat is extracted with a mixture of acetonitrile/water, centrifuged, and chilled. The acetonitrile phase is diluted with 2% NaCl and the synthetic pyrethrins are partitioned into hexane. A portion of the hexane extract is further purified using alumina column chromatography. The column eluate is concentrated, adjusted to 6% acetonitrile in phosphate-buffered saline, and analyzed with a competition enzyme-linked immunosorbent assay (c-ELISA).



#### 2. Applicability

This method is applicable to bovine, porcine, and poultry fat. Synthetic pyrethrins are used for their insecticidal action. The monoclonal antibody used in this assay will detect the following pyrethroid insecticides: permethrin, cypermethrin, deltamethrin, phenothrin, and fenpropathrin. Flucythrinate and fenvalerate are only weakly detected, and tetramethrin is not recognized by the antibody. Refer to section J.7 for sensitivities to various pesticides. Identification and quantitation of the individual pyrethroids detected is not possible, because ELISA response is integrated over the spectrum of all detectable compounds.

# SYNTHETIC PYRETHRINS/C-ELISA

# **DETERMINATIVE METHOD**

## **B. EQUIPMENT**

1. Apparatus	a.	Centrifuge (Sorval T6000B with adapter for 50 mL tubes).
	b.	Centrifuge tubes, 50 mL polyproylene disposable, with cap (Beckton- Dickinson Falcon 2098).
	С.	Water bath, heated circulating (Freas model 260).
	d.	125 mL glass separatory funnel with stopper.
	е.	Centrifuge tube, glass, 15 mL graduated, with glass stopper.
	f.	Pasteur pipet, disposable, 5.25" (Scientific Products, Cat. P5202-1).
	g.	Pipetter, 50-200 $\mu$ L adjustable, 8-channel, with tips (Titertek digital multichannel pipette, Flow Laboratories).
	h.	Pipetters, 2-20, 10-100, and 100-1000 $\mu$ L adjustable single-channel, with tips (Gibson Pipetman P-20, P-100, P-1000).
	i.	Immulon II 96-well microtiter plates (Dynatech Laboratories).
	j.	Sealing tape, 8.26 $\times$ 13.34 cm, for microtiter plates (Dynatech Laboratories, Cat. 001-010-3501).
	k.	Skatron 96-well microtiter plate washer (Skatron, Inc.).
	1.	Incubator, mechanical convection (Precision Scientific Model 4EM).
	m.	Soxhlet extraction apparatus (Kontes 585000-023).
	n.	Extraction thimble, cellulose, 43 $\times$ 123 mm (Whatman 2800432).
	о.	Filter paper, Whatman #1, 24 cm.
	NC	DTE: Equivalent apparatus may be substituted for those specified.
2. Instrumentation	a.	Bio-Tek Model 310-96 well-plate reader with software (Bio-Tek Instruments, Inc.).

b. IBM PC/AT or compatible computer.

## PYRE-2

#### C. REAGENTS AND SOLUTIONS

- a. Acetonitrile (ACN).
- b. Hexane.
- c. Benzene.
- d. Dichloromethane (DCM).
- e. Polyoxyethylene-sorbitan (Tween 20, Sigma Chemical Co.).
- f. Sodium chloride, reagent grade.
- g. Potassium phosphate, monobasic, reagent grade.
- h. Potassium phosphate, dibasic, reagent grade.
- i. Sodium azide (Sigma Chemical Co.).
- j. Magnesium chloride hexahydrate.
- k. Diethanolamine, 98% (Sigma Chemical Co.).
- I. p-Nitrophenyl phosphate, 5 mg tablets (Sigma Chemical Co.). Store in a freezer at -10° C.
- m. Alumina oxide (Bio-Rad acid alumina AG-4 100-200 mesh). Extract with dichloromethane for 24 hours in a soxhlet apparatus. Dry at 130° C for 24 hours; then store at 110° C.
- n. Ovalbumin, crude powder, grade II (Sigma Chemical Co.).
- Anti-mouse IgG (whole molecule) alkaline phosphatase conjugated, affinity isolated antigen, specific antibody developed in goat (Sigma Chemical Co.). Store at 4° C.
- p. Py-1 antibody, 1 mg/mL. Monoclonal anti-pyrethroid antibody is referred to as Py-1 (Stanker, L. H., et al., *J. Agr. Food Chem.* 37:834-839, 1989). Py-1 is produced by the corresponding hybridoma cell line as an ascites fluid. The antibody is purified and stored at  $-20^{\circ}$  C as a stock solution at a concentration of 1 mg/mL in small aliquots. Once an aliquot is thawed, it should be stored at 4° C. When used in the competition ELISA described in this method, using permethrin as a competitor, an I<sub>so</sub> of 1.55 ng  $\pm$  0.6 ng should be observed. (For source of supply, contact Chemistry Division, FSIS.)
- q. Coating antigen stock solution, 5 mg/mL: 3-phenoxyacetic acid conjugated to bovine serum albumin (3-pba-BSA) in PBS-7. The quality of each lot of conjugate must be evaluated in a c-ELISA using permethrin as a competitor and the monoclonal antibody Py-1. An I₅₀ of 1.55 ng ± 0.6 ng should be observed. Store at 20° C in small aliquots. (For source of supply, contact Chemistry Division, FSIS.)

NOTE: All solvents should be certified for use in pesticide analysis.

## SYNTHETIC PYRETHRINS/C-ELISA

PYRE-3

1. Reagent List

# FSIS

# **DETERMINATIVE METHOD**

## C. REAGENTS AND SOLUTIONS (Continued)

2. Solutions List	а.	Tween 20, 1% solution. Dilute 1 mL Tween 20 to 100 mL with distilled water.
	b.	Tween 20, 0.05% wash solution. Dilute 1 mL of Tween 20 to 2 L with distilled water.
	C.	Phosphate buffered saline, pH 7.0 (PBS-7). Combine 9.65 g Na <sub>2</sub> HPO <sub>4</sub> , 1.9 g NaH <sub>2</sub> PO <sub>4</sub> , and 42.5 g NaCl. Add 9 L of distilled water and ajdust the pH to 7.0 using 1N HCl or NaOH. Dilute to 10 L with distilled water.
	d.	Assay buffer. Dilute 1 mL of 1% Tween 20 solution to 1 L with PBS-7 and mix thoroughly. Store at 4° C.
	e.	Phosphate buffered saline, pH 9.0 (PBS-9). Add 61.4 g NaCl and 18.76 g Na <sub>2</sub> HPO <sub>4</sub> to 6 L of distilled water, and adjust the pH to 9.0 using 1N NaOH. Adjust volume to 7 L. Store at 4° C.
	f.	Diethanolamine buffer, 10%. Add 97 mL of diethanolamine, 0.2 g NaN <sub>3</sub> , and 0.1 g MgCl <sub>2</sub> •6H <sub>2</sub> O to 800 mL distilled water. Adjust pH to 9.8 with 6N HCl and adjust final volume to 1 L. The buffer is filtered (Whatman #1) and stored at 4° C.
	g.	pba-BSA, 5 $\mu$ g/mL solution. Dilute 100 $\mu$ L of the 5 mg/mL pba-BSA stock solution to 100 mL with distilled water. This provides enough solution to coat ten 96-well plates. Prepare fresh prior to coating plates.
	h.	Ovalbumin, 3% in PBS-9. Add 4.5 g of ovalbumin powder to a 200-250 mL flask or bottle containing 150 mL of PBS-9 buffer at 30-40° C, cap, and mix until dissolved. Filter through glass wool and collect in a stoppered bottle. Store at 4° C.
	i.	Py-1 antibody, 1:5000 dilution. Dilute 3 $\mu L$ Py-1 antibody (1 mg/mL) to 15 mL with assay buffer. Prepare immediately before use.
	j.	Antimouse IgG, 1:500 dilution in assay buffer. Dilute 30 $\mu L$ of antimouse IgG, 5 mg/mL, to 15 mL with assay buffer. Prepare immediately before use.
	k.	p-Nitrophenyl phosphate solution, 1 mg/mL. Dissolve three 5 mg tablets of p-nitrophenyl phosphate in 15 mL 10% diethanolamine buffer. Prepare immediately before use.
	I.	ELISA control solution. 6% ACN in assay buffer. Dilute 6 mL acetonitrile to 100 mL with assay buffer.
	m.	Extraction solvent, 85% acetonitrile/water. Add 850 mL acetonitrile to a 1 L volumetric flask or mixing cylinder and dilute to volume with distilled water.
	n.	2% NaCl solution. Add $20$ g NaCl to a 1 L volumetric flask or mixing cylinder and dilute to volume with distilled water.
	0.	6% acetonitrile/assay buffer. Add 6.0 mL of acetonitrile to a 100 mL volumetric flask and dilute to volume with assay buffer.

#### C. REAGENTS AND SOLUTIONS (Continued)

3. Preparation of Microtiter Plates	a.	Using a multichannel pipet, add 100 $\mu$ L of a 5 $\mu$ g/mL solution of pba-BSA in distilled water to each well of a 96-well microtiter plate.
	h	Place the plate uncovered in an incubator of 27% C and allow the

- b. Place the plate, uncovered, in an incubator at 37° C, and allow the pba-BSA solution to evaporate overnight. Make sure that all traces of moisture are gone before removing plate.
- c. Store "coated" plates in a sealed container at 4° C, and use within 4 weeks.

#### D. STANDARDS

1. Source	Permethrin (mixed isomers) and other synthetic pyrethroids may be obtained from:
	Director Environmental Monitoring Support Laboratory Office of Research and Development Cincinnati, OH 45268
2. Preparation of Standards	a. Permethrin stock standard, 0.5 mg/mL permethrin. Accurately weight out 25 mg of the permethrin standard (mixed isomers) into a 50 mL beaker and dissolve in ca. 30 mL acetonitrile. Transfer quantitatively to a 50 mL volumetric flask and dilute to volume with acetonitrile.
	b. Permethrin intermediate standard, 10 μg/mL. Pipet 1.0 mL of permethrin stock standard into a 50 mL volumetric flask and dilute to volume with acetonitrile.
	c. Permethrin fortification standard, 2.5 $\mu$ g/mL. Pipet 5 mL of 10 $\mu$ g/mL permethrin standard into a 25 mL volumetric flask. Pipet 15 mL acetonitrile into the flask (20 mL final volume) and mix.
	d. Permethrin ELISA standard, 0.3 μg/mL. Prepare a 0.30 μg/mL standard to be used in constructing a standard curve by adding 60 μL of the 10 μg/mL permethrin standard into a 15 mL centrifuge tube and diluting to 2 mL with assay buffer.
3. Storage Conditions	Store all standards at 4° C.
4. Shelf Life Stability	a. Stock, intermediate, and fortification standards are stable for at least 6 months.
	b. Permethrin ELISA standard must be prepared fresh daily.

-

# **DETERMINATIVE METHOD**

## E. EXTRACTION PROCEDURE

1. Sample Preparation	Pas gla the to (	ss chilled sample through a food chopper twice. Plug a powder funnel with ss wool. Put approximately 100 g of the comminuted fat into the funnel, place funnel in a beaker, and render at 100-110° C overnight or until the fat ceases drop.
2. Sample Extraction	a.	Add 2.5 g of rendered fat into a 50 mL centrifuge tube. Add 25 mL 85% acetonitrile/water extraction solvent, cap securely, and place tube into a 55° C water bath for 5 min.
	b.	Shake sample vigorously for 3 min. A mechanical shaker may be used for this operation if it has a cycle speed greater than 200 strokes/min.
	c.	Immediately centrifuge sample for 2 min at 3000 rpm. Chill sample until fat has solidified. (Do not chill to point of phase separation or freezing of water phase.)
	d.	Remove 10 mL of the liquid phase and place in a 125 mL separatory funnel. No suspended particles of fat should be transferred.
	e.	Add 50 mL 2% NaCl solution and swirl. Add 15.0 mL hexane, stopper, and shake for 2 min.
	f.	Carefully drain and discard the aqueous fraction.
	g.	Add 10 mL of distilled water to the separatory funnel and shake gently for 30 sec. Allow phases to separate completely.
	h.	Drain and discard the 10 mL wash. A small amount of hexane may be drained to displace residual water in the stopcock.
	i.	Drain the hexane phase into a 15 mL graduated centrifuge tube. Adjust volume to 15 mL with hexane.
	j.	Transfer the hexane eluate to a 1 cm $\times$ 0.5 cm id column of alumina topped with 2 cm sodium sulfate. Avoid transferring any residual water that may be in the tip of the centrifuge tube. If less sensitivity is required, a suitable aliquot of the eluate may be diluted to 15 mL and applied to the column.
	k.	Wash the column with two 5 mL portions of hexane. Discard the washes.
	١.	Elute permethrin from the column into a 15 mL centrifuge tube using 5 mL benzene.
	m.	Evaporate the organic solvent to dryness under a stream of nitrogen in a 40° C water bath. Dissolve the residue in 60 $\mu$ L of acetonitrile, then dilute to 1.0 mL with assay buffer.
	n.	Analyze sample extracts for permethrin concentration using the competition ELISA described in section F.

## **DETERMINATIVE METHOD**

#### E. EXTRACTION PROCEDURE (Continued)



#### F. ANALYTICAL QUANTITATION

1. Competition ELISA	NOTE: All solutions used should be brought to room temperature before being
·	added to microtiter plates.

- a. Add 400  $\mu$ L of 3% ovalbumin solution to each well of an antigen-coated microtiter plate. Incubate the plate at 25-37° C for 1 hour. Wash the plate 3 times with 0.05% Tween 20/water solution. Remove residual wash solution by inverting the plate and rapping repeatedly onto an absorbent surface (e.g., paper towels).
- b. Refer to the plate map in section F.2 below. Add 100  $\mu$ L of 6% acetonitrile in assay buffer to columns 1, 2, and 12, and to all wells in row A except A3. Add 200  $\mu$ L of 0.3  $\mu$ g/mL standard solution to well A3.
- c. Prepare a standard curve by making seven serial 1:2 dilutions. Transfer 100  $\mu$ L of standard from well A3 to well A4, and mix by repeated pipetting. Repeat the procedure for wells A4-A5, A5-A6, A6-A8 (A7 is a control), A8-A9, A9-A10, and A10-A11. Remove 100  $\mu$ L from A11 and discard. (Standard in well A3 is equivalent to 300 ppb permethrin in the sample unless an aliquot is taken in section E.2, step j.
- d. All samples are analyzed in triplicate. Using the plate map in F.2 as a template, add 100  $\mu$ L aliquots of each sample to each of the three specified wells. After all samples have been applied, any empty wells remaining can be filled with 100  $\mu$ L of 6% acetonitrile in assay buffer. These will serve as additional controls.
- e. Add 100  $\mu$ L of a 1/5000 dilution of the 1 mg/mL stock solution of Py-1 in assay buffer to all wells except those in column 1. Column 1 is the no primary antibody blank.
- f. Seal the plate using a plate sealer or other airtight wrap and incubate for 1 hour at 37° C.
- g. Wash the plate 7 times with 0.05% Tween 20 solution. Remove residual wash solution as in step a above.
- h. Pipet 50  $\mu$ L of alkaline phosphatase conjugated goat-antimouse immunoglobulin diluted 1/500 in assay buffer into all wells of the plate.
- i. Seal and incubate the plate at 37° C for 1 hour; then wash as in step g.
- j. Add 100  $\mu$ L of a 1 mg/mL solution of p-nitrophenyl phosphate (freshly prepared in 10% diethanolamine buffer) to all wells. Allow color to develop at room temperature for 30 minutes, or until the absorbance of the controls is in the range 1.0-2.0; then record the absorbances of all wells in dual wavelength mode at 405 nm (540 nm reference) using a plate reader.

## SYNTHETIC PYRETHRINS/C-ELISA

## **DETERMINATIVE METHOD**

## F. ANALYTICAL QUANTITATION (Continued)

#### 2. ELISA Plate Distribution Map

		-	-		_	-		-				
	1	2	3	_ 4	5	6	7	8	9	10	11	12
A	В	С	S	S/2	S/4	S/8	С	S/16	S/32	S/64	S/128	С
В	В	С	1	2	3	4	5	6	7	8	9	С
С	В	С	10	11	12	13	14	15	16	17	18	С
D	В	С	19	20	21	1	2	3	4	5	6	С
E	В	С	7	8	9	10	11	12	13	14	15	С
F	В	С	16	17	18	19	20	21	1	2	3	С
G	В	С	4	5	6	7	8	9	10	11	12	С
H	В	С	13	14	15	16	17	18	19	20	21	С

- B = Blanks (no antibody; 6% acetonitrile/assay buffer)
- C = Controls (6% acetonitrile/assay buffer)
- S = Standard (30 ng/well)
- 1-21 = Samples applied to plate (each sample is applied three times)

#### F. ANALYTICAL QUANTITATION (Continued)

3. Competition ELISA Flow Chart



## SYNTHETIC PYRETHRINS/C-ELISA

PYRE-11

**FSIS** 

## DETERMINATIVE METHOD

#### G. CALCULATIONS

#### 1. Procedure

NOTE: Since the ELISA antibody integrates its response over the entire spectrum of compounds that it recognizes, it is not possible to determine the identity or actual concentration of the pyrethroids present in a positive sample. All positives are quantitated as ppb permethrin.

a. Each row on the plate is treated as a separate experiment. Determine an average control value for each row by averaging all acceptable control values in the row. Any individual value that appears to be an outlier may be tested and discarded. Once average values have been determined, the percent inhibition for all wells in columns 3-6 and 8-11 can be calculated using the formula:

Percent inhibition = [Abs. (control) - Abs. (sample)] Abs. (control) - Abs. (blank) X 100%

- b. Average the three separate values of percent inhibition for each sample. Any individual value that appears to be an outlier may be tested and discarded.
- c. Using semi-log paper, plot concentrations of the permethrin standards in ng/well (X axis) versus their percent inhibition values. Connect all points to form a standard curve. Concentrations (as permethrin) of all samples considered positive (see NOTE below) can be determined graphically by constructing a drop line from the sample's percent inhibition value on the standard curve to the X axis.

Convert concentrations to ppb permethrin by dividing the permethrin concentrations in ng/well by the sample weight in grams/well. If no dilutions are made in section E.2, step j, this value is 0.1 g. If dilutions were made, this value may be calculated using the formula:

g sample/well = 2.667(W)(A)/1000V, where

- W = sample weight in grams A = mL extract from E.2.i applied to alumina column
- V = final volume of extract in mL

NOTE: It is not uncommon for blank samples to exhibit high background readings. Percent inhibition readings as high as 15-20% have been observed with some blank fat samples. Samples giving inhibition values at or below this level range are likely to show a high incidence of false positives. Reliable detection of probable positive samples requires > 25% inhibition of response.

## G. CALCULATIONS (Continued)

# 2. Standard Curves



**Standard Curve for Permethrin** 

Permethrin (PPB)

# SYNTHETIC PYRETHRINS/C-ELISA

**PYRE-13** 

**FSIS** 

## **DETERMINATIVE METHOD**

## G. CALCULATIONS (Continued)



Fortified Standard Curve in Poultry Fat

Permethrin (PPB)

## G. CALCULATIONS (Continued)



Fortified Standard Curve in Swine Fat

Permethrin (PPM)

# SYNTHETIC PYRETHRINS/C-ELISA

**PYRE-15** 

**FSIS** 

# **DETERMINATIVE METHOD**

## G. CALCULATIONS (Continued)



**Standard Curves for Synthetic Pyrethroids** 

**Pyrethrins (PPB)** 

**SYNTHETIC PYRETHRINS/C-ELISA** 

----

# **DETERMINATIVE METHOD**

## H. HAZARD ANALYSIS

1. Method Title	Immunoassay Determina	tion of Synthetic Pyrethrins	by C-ELISA		
2. Required Protective Equipment	Safety glasses, plastic gloves, lab coat.				
3. Procedure Steps		Hazards	Recommended Safe Procedures		
	C. Reagents				
	Benzene Acetonitrile Hexane	These solvents are all flammable and can produce toxic effects through exposure to the skin, eyes, respiratory system, or ingestion. Benzene is a carcinogen.	All procedures involving solvents should be carried out in an efficient fume hood and away from any source of ignition. Protective gear should be worn at all steps where the possibility of solvent spillage exists.		
	D. Standards	Concentrated standards may be toxic. Exposure paths are the same as described above for reagents.	Prepare standards in a fume hood or glove box. Use protective gear.		
4. Disposal Procedures	Reagent mixtures used in section E	(see above)	Hold in designated cans awaiting disposal by the contractor or in-house specialist.		
_	Alumina columns (used)	(see above)	Allow columns to dry in a fume hood. Dispose of packing by pouring into a plastic bag and discarding.		

# **DETERMINATIVE METHOD**

## J. QUALITY ASSURANCE PLAN

1. Performance Standards	Compound	Acceptable Recovery (%)	Analytical Range (ppb)	Repeatability %CV	
	Permethrin	50-100	50-200	< 30	
2. Critical Control Points and	Record	1	Acceptable	Control	
Specifications	Reagents, solvents (C	.1) N	Must be free of interfering substances		
	Sample weight (E.2.a)	) 2	2.50 ± 0.05 g		
	Column elution rate (E	E.2.j-l) L	Less than 1.0 mL/minsteps	n for all elution	
	Incubation temperatur	re (F.1.g,j)	37° ± 3° C		
	Incubation conditions	(F.1.g,j) f i	Plates must be tightly sealed prior to incubation steps		
3. Readiness To Perform	a. Familiarization. i. Phase I: Star 200 ppb with ii. Phase II: Ber	ndards—Analyze 1 three replicates	permethrin standard for each compound	s at 50, 100, and ude at least three	
	levels with the permethrin st tissue blank.	andards. Each s	ach at 50, 100, and et should include a r	eagent blank and	
	NOTE: Phase	e I and Phase II	may be performed co	oncurrently.	
	iii. Phase III: Ch each species range 50-200	analyst accreditation ding blanks and tissu	-14 samples for ues fortified in the		
	(a) Samples	s submitted by su	upervisor or LSO.		
	(b) Report (	data to Chemistry	y Division, QSB.		
	Notification from C	hemistry Division	required to commenc	e official analysis.	
	b. Acceptability crite	ria.			
	i. No false pos	itives.			
	ii. No recovery	missed if preser	nt at or above 50 pp	b (fortified level).	
	iii. Refer to sect	tion J.1 above.			

.....

# DETERMINATIVE METHOD

J. QUALITY ASSURANCE	E PL/	AN (Continued)
4. Intralaboratory	a.	System, minimum contents.
Check Samples		i. Frequency: At least one check sample biweekly per analyst.
		<li>Blind check samples or random duplicates chosen by the supervisor or LSO. Records to be maintained by the analyst and reviewed by the supervisor and LSO:</li>
		(a) All duplicate findings.
		(b) All recovery values.
		(c) Running average, standard deviation, and CV for all recoveries.
		(d) CUSUM charts.
	b.	Acceptability criteria.
		If unacceptable values are obtained, then:
		i. Stop all official analyses for that analyst.
		ii. Investigate and identify probable cause.
		iii. Take corrective action.
		iv. Repeat Phase II of section J.3 above if cause was analyst-related.
5. Sample Acceptability	a.	Matrix: Fat.
and Stadility	b.	Sample receipt size: Sufficient tissue for all quantitative and confirmational analyses and sample reserve portion (approximately 100 g).
	c.	Sample receipt condition: Chilled to frozen.
	d.	Sample storage:
		i. Time: 3 months.
		ii. Condition: Store at -4° C or below until analyzed.
6. Sample Set	Ea	ch set must contain a tissue blank, a fortified blank at 50-200 ppb (100 ppb commended), and the samples to be analyzed.
7. Sensitivity	a.	Lowest detectable limit (LDL), ppb: permethrin (mixed), 25; others, see 7.c.
	b.	Lowest reliable quantitation (LRQ), ppb: permethrin (mixed), 50; others, see $7.c.$

# SYNTHETIC PYRETHRINS/C-ELISA

# PYRE

July, 1991

## **DETERMINATIVE METHOD**

### J. QUALITY ASSURANCE PLAN (Continued)

c.	Py-1 antibody sensitivity	for selected pyrethroi	ds.
	Compound	I₅₀ Value (ng)	Method LDL (ppm)
	Permethrin (mixed)	1.5	0.025
	Phenothrin	1.5	0.025*
	Deltamethrin	15	0.25*
	Fenpropathrin	17	0.28*
	Cypermethrin	30	0.50*
	Flucythrinate	120	2.0*
	Fenvalerate	350	5.8*
	Tetramethrin	> 600	> 10*