

University of Manchester - Chondrocyte Protocol (Susan Kimber Lab)

Title	University of Manchester - Chondrocyte protocol
Date Submitted	Jan 2012
Submitted by -	Sunita DSouza
Adapted from -	Directed differentiation of human embryonic stem cells toward chondrocytes Rachel A Oldershaw , Melissa A Baxter , Emma T Lowe , Nicola Bates , Lisa M Grady , Francesca Soncin , Daniel R Brison , Timothy E Hardingham & Susan J Kimber Nature Biotechnology 28, 1187–1194 (2010) Published online, 22 October 2010
Contributors -	Rachel A Oldershaw , Melissa A Baxter , Emma T Lowe , Nicola Bates , Lisa M Grady , Francesca Soncin , Daniel R Brison , Timothy E Hardingham & Susan J Kimber
University of Manchester	University of Manchester, UK

This next section will be on the side and is only relevant for website programmers

Score	=Likes/(Likes+Dislikes)+Likes/Downloads or some other algorithm... Suggestions??? I can also come up with something better in due time.
Downloads	#
“Likes”	#
“Dislikes”	#
Status	[Validated (>=3, In Progress (=1-2), Not Validated (none)]
Validated Core 1	Name / Institute
Validated Core 2	Name / Institute
Validated Core 3	Name / Institute


Table of Contents

1. Introduction
2. FlowChart
3. Materials
 - a) Reagents
 - b) Media Preparation
4. Protocol
 - a) Feeder – free culture
 - b) Directed Differentiation Protocol
5. Acknowledgements
6. Contributors


INTRODUCTION

We have developed for hESC a new 3-Stage directed differentiation protocol (DDP) to generate chondrocytes, the specialized cells that form cartilage tissue. The protocol is segmented into stages that mimic the developmental processes that occur in cell lineage specification during embryogenesis. Pluripotent hESC, are established as feeder-free cultures⁴ prior to DDP being carried out in monolayer culture on a sequence of matrix protein-coated surfaces with a chemically defined medium sequentially supplemented with growth factors to direct differentiation through developmental intermediate cell populations of primitive streak/mesendoderm (Stage 1), mesoderm (Stage 2) and then chondrocytes (Stage 3) (Figure 1). This unique protocol is highly efficient, scalable and completely chemically-defined, thus making it appropriate for translation towards clinical applications of chondrocytes, as well as providing a defined system for characterizing the molecular mechanisms regulating hESC differentiation.

FLOWCHART:

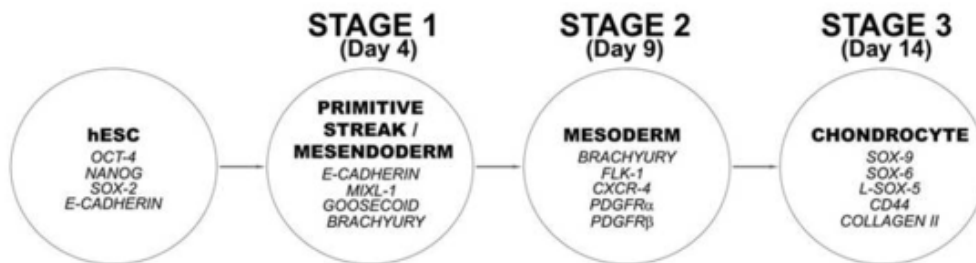

FIGURES

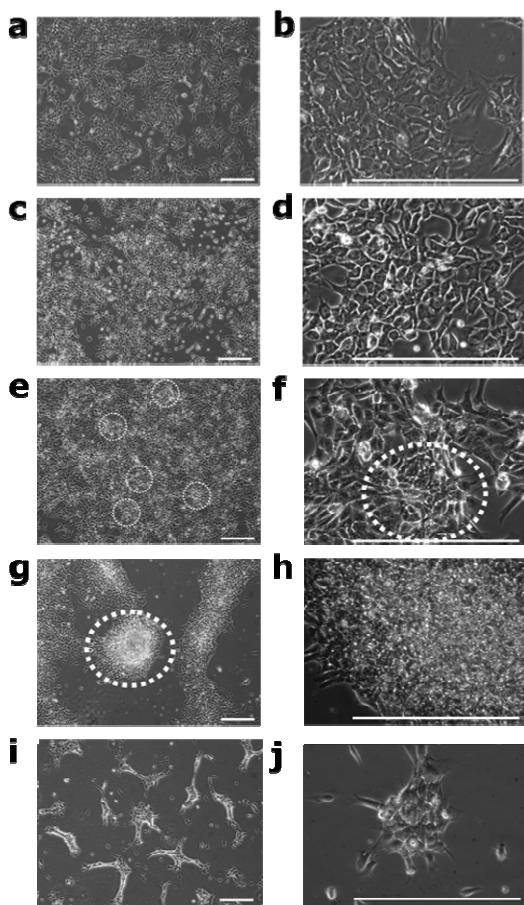

Cell culture plastics:

	Company	Cat #
35mm Culture Dishes BD BioCoat™ Poly-L-Lysine	BD Biosciences	354518
6 Well Costar® Clear TC-Treated Multiple Well Plates	Corning	3516
5ml Costar® Stripette® polystyrene serological pipettes	Corning	4487
10ml Costar® Stripette® polystyrene serological pipettes	Corning	4488
25ml Costar® Stripette® polystyrene serological pipettes	Corning	4489
20µl Bevelled Filter Tip	Starlab	S1120-1810
200µl Graduated Filter Tips	Starlab	S1120-8810
1000µl Filter Tip	Starlab	S1126-7810
1.5ml Microcentrifuge Tube	Starlab	E1415-1500
15ml polypropylene centrifuge tubes	Corning	CLS430790
50ml polypropylene centrifuge tubes	Corning	CLS430790
0.22µm syringe-driven Millex-GP Filter Unit	Millipore	SLGP033RB
0.22µm vacuum-driven Steritop-GP Filter Unit	Millipore	SCGPT01RE
Stericup Receiver Flask, 45 mm thread	Millipore	SC00B05RE
20ml syringe	VWR	613-3921
Knockout™ DMEM	Invitrogen	10829-018
Knockout™ Serum Replacement	Invitrogen	10828-028
DMEM:F12 high glucose with 2mM L-glutamine and 15mM HEPES buffer	Invitrogen	11330-032
DMEM high glucose	Invitrogen	31053-028
Fetal Bovine Serum	Invitrogen	10082-147
Insulin Transferrin Selenium (ITS) Supplement (100X)	Invitrogen	51300-044
MEM Non-Essential Amino Acids (NEAA) 10mM (100X)	Invitrogen	11140-035
L-Glutamine 200mM (100X) (To prevent freeze-thaw cycles the L-Glutamine is thawed, aliquoted in 5ml volumes and stored at -20 °C)	Invitrogen	25030-024
Penicillin-Streptomycin (100X) contains 5,000 units of penicillin (base) and 5,000 µg of streptomycin (base)/ml utilizing penicillin G (sodium salt) and streptomycin sulfate in 0.85% saline (To prevent freeze-thaw cycles the Penicillin-Streptomycin is thawed, aliquoted in 5ml volumes and stored at -20 °C)	Invitrogen	
β-Mercaptoethanol 55mM in DPBS	Invitrogen	31350010
N2 Liquid Supplement (100X)	Invitrogen	17502-048
B27 Liquid Supplement (50X)	Invitrogen	17504-044
Dulbecco's Phosphate Buffered Saline (DPBS) (1X)	Invitrogen	14040-091
Trypsin:EDTA (To prevent freeze-thaw cycles the Trypsin:EDTA is thawed, aliquoted in 10ml volumes and stored at -20 °C)		L11-660
Distilled H2O (dH2O) is sterilized by autoclaving	-	-
Dimethyl Sulfoxide	Sigma	D2650
Porcine gelatin	Sigma	G1890
Bovine Serum Albumin (BSA)	Sigma	A4161
Human Plasma Fibronectin	Millipore	FC010
Human recombinant FGF2	Peptotech	AA10-155

Human recombinant Activin-A	Peprotech	120-14
Human recombinant Neurotrophin-4	Peprotech	450-04
Human recombinant BMP4	Peprotech	120-05
Human recombinant GDF5	Peprotech	BMP14/CDMP1
Human recombinant Follistatin 300	Peprotech	F1175-25UG
Mouse recombinant Wnt3a	Peprotech	1324-WN

Tip: Due to the decrease in the biological activity of supplementary growth factors we store all culture media at 4°C and discard 14 days after preparation. In order to prevent waste from discarding expired media we frequently prepare 500ml volumes of hESc Feeder-Free Medium and Directed Differentiation Basal Medium without the growth factors and store 40ml aliquots at -20°C. The required growth factor combinations are then added to thawed aliquots. We have not detected any adverse effects in cell cultures when using media that has been stored at -20°C for up to 2 months from preparation.

Growth factors:

Growth factors are purchased as lyophilized powders which are reconstituted to a stock concentration with 0.1% (vol/vol) BSA in DPBS. To prepare 0.1% (vol/vol) BSA add 0.01g of BSA to 10ml DPBS, vortex well and sterilize through a 0.22µm syringe driven filter.

Human recombinant FGF2 (AA10-155) (Biosource™, Invitrogen cat. no. PHG0021) – Dissolve 1mg of lyophilized powder in 10ml of 0.1% (vol/vol) BSA in DPBS to make a stock concentration of 100µg/ml (wt/vol). Aliquot the stock growth factor solution in 50µl volumes.

Human recombinant Activin-A (Peprotech cat. no. 120-14) – dissolve 5µg of lyophilized powder in 500µl of 0.1% (vol/vol) BSA in DPBS to make a stock concentration of 10µg/ml (wt/vol). Aliquot the stock growth factor solution in 50µl volumes.

Human recombinant Neurotrophin-4 (NT4) (Peprotech cat. no. 450-04) – Dissolve 10µg of lyophilized powder in 1ml of 0.1% (vol/vol) BSA in DPBS to make a stock concentration of 10µg/ml (wt/vol). Aliquot the stock growth factor solution in 20µl volumes.

Human recombinant BMP4 (Peprotech cat. no. 120-05) – Dissolve 5µg of lyophilized powder in 500µl of 0.1% (vol/vol) BSA in DPBS to make a stock concentration of 10µg/ml (wt/vol). Aliquot the stock growth factor solution in 50µl volumes.

Human recombinant GDF5 (BMP14/CDMP1) (Peprotech cat. no. 120-01) – Dissolve 50µg of lyophilized powder in 5ml of 0.1% (vol/vol) BSA in DPBS to make a stock concentration of 10µg/ml (wt/vol). Aliquot the stock growth factor solution in 50µl volumes.

Human recombinant Follistatin 300 (Sigma cat. no. F1175-25UG) – Dissolve 25µg of lyophilized powder in 2.5ml of 0.1% (vol/vol) BSA in DPBS to make a stock concentration of 10µg/ml (wt/vol). Aliquot the stock growth factor solution in 50µl volumes.

Mouse recombinant Wnt3a (R&D Systems cat. no. 1324-WN) – Dissolve 2µg of lyophilized powder is dissolved in 200µl of 0.1% (vol/vol) BSA in DPBS to make a stock concentration of 10µg/ml (wt/vol). Aliquot the stock growth factor solution in 25µl volumes.

Critical: Growth factor polypeptides are highly unstable and biological activity decreases rapidly. Because of this lyophilized growth factors are assumed to be sterile on opening and reconstitution is carried out in a microbiological safety cabinet using aseptic technique. Growth factor solutions are never passed through 0.22µm filters and cell culture media are sterilized prior to the addition of growth factors. Once reconstituted the growth factor aliquots are transferred on ice for storage at -80°C. Thawed aliquots are never re-frozen but are stored at 4°C and discarded 14 days post-thaw.

Cryopreservation medium: Add a 1ml volume of Dimethyl Sulfoxide (Sigma cat. no. D2650) to 9ml volume of FBS. Pipette up and down to mix well.

Porcine gelatin (Sigma cat. no. G1890) – A 1% stock gelatin solution is prepared by adding 4g of lyophilized gelatin powder to 400ml dH₂O. The solution is autoclaved to sterilise, aliquotted in 20ml volumes and stored at -20°C. Working gelatin solutions of 0.1% (vol/vol) are prepared in sterile T-75 cell culture flasks by adding 20ml of thawed 1% (vol/vol) gelatin to 180ml of sterile dH₂O. The 0.1% (vol/vol) gelatin solution is stored at 4°C and discarded after 28 days

Bovine Serum Albumin (BSA) (Sigma cat. no. A4161)
 Human Plasma Fibronectin (Millipore cat. no.FC010) – Fibronectin is supplied as a liquid at a concentration of 1mg/ml (wt/vol) to be stored at 4°C. Working solutions of 50µg/ml (wt/vol) are prepared by diluting 1ml of stock fibronectin in 19ml of DPBS and stored at 4°C.

Table 1: Formulation of MEF Medium

MEF Medium (500ml)	
DMEM	440ml
FBS	50ml
L-Glutamine	5ml
Penicillin-Streptomycin	5ml
Pass the medium through a 0.22µm filter and store at 4°C for up to 14 days	

Table 2: Formulation of hESc Medium

hESc Medium (200 ml)	
Knockout™ DMEM	154ml
Knockout™ Serum Replacement	40ml
ITS	2ml
NEAA	2ml
L-Glutamine	2ml
β-Mercaptoethanol	367µl
Pass the medium through a 0.22µm filter and add 20µl of 100µg/ml (wt/vol) stock FGF2. Store at 4°C for up to 14 days	

hESc culture on Inactivated Mouse Embryonic Fibroblast (iMEF) Feeder Cell Layer:
 1. Transfer 1ml of 0.1% (vol/vol) gelatin to 1×35mm culture dish and incubate in a cell culture incubator at 37°C for 1 hour.

2. Remove one cryovial of iMEF cells from the liquid nitrogen cell storage dewar and place on ice for transfer to the cell culture suite.
3. Hold the cryovial by the lid and submerge three quarters of the cryovial in the 37°C water bath for 2-3 minutes to thaw the cell suspension. **Tip: Swirl the cryovial in the water to facilitate heat transfer. Thaw the iMEF cells until approximately 10% of ice remains, the iMEF cells will continue to thaw during transfer to the microbiological safety cabinet.**
4. Draw up 4ml of MEF medium in a 5ml serological pipette and then the 1ml volume of iMEF cell suspension from the cryovial. Transfer the 5ml volume to a sterile 15ml polystyrene centrifuge tube.
5. Add 90µl of the iMEF cell suspension and 10µl of trypan blue solution to a 1.5ml microfuge tube and mix well by pipetting up and down. Transfer 10µl to each chamber of haemocytometer and perform a viable cell count. MEF cells are seeded at a cell density of 3×10^4 cells/cm² which for a 35mm culture dish is 2.92×10^5 cells; calculate the volume of iMEF cell suspension required: 2.92×10^5 cells / viable cells/ml = Xml of MEF cell suspension. Transfer the calculated volume to a sterile 15ml polystyrene centrifuge tube and centrifuge at 720 x g for 3 minutes at room temperature. Remove the supernatant and resuspend the cell pellet in 2ml of MEF medium (Table 1).
6. Take the 35mm culture dish with gelatin from the incubator, remove the gelatin and add the 2ml iMEF cell suspension. Transfer to the cell culture incubator.
7. Twenty-four hours post-seeding check the iMEF cells under the microscope to ensure that they have adhered to the cell culture surface and formed a homogeneous culture of approximately 80-85% confluence.
8. Thaw hESc and seed onto the iMEF layer. Remove a vial of hESc from liquid nitrogen storage and thaw as for the iMEF cells in steps 3-5 but substitute hESc medium (Table 2) for MEF medium. Centrifuge the hESc at 720 x g for 3 minutes at room temperature and resuspend the cell pellet in 2ml of hESc Medium.
9. Remove the iMEF medium from the iMEF cell layer and add the hESc suspension. Transfer the cell culture dish to the cell culture incubator.
10. Change the hESc medium every 2 days and check the hESc for growth under the microscope. Colonies of hESc can be observed growing between iMEF cells after approximately 2-4 days.
11. Passage the hESc at a ratio of 1:6 when the culture has grown to approximately 90-95% confluence. Prepare 6 wells of a 6 well culture dish with a iMEF feeder cell layer as described in steps 1-7.
12. Pre-warm a 10ml aliquot of Trypsin:EDTA and hESc medium by placing in the 37°C water bath for 30 minutes.
13. Remove and discard the hESc medium from the hESc culture and add 1ml of Trypsin:EDTA. Incubate in the cell culture incubator at 37°C and check under the microscope every 2-3 minutes for cell dissociation. iMEF cells are seen to round up whereas hESc can be seen to separate from each other. **Critical: Cell:cell contact is crucial in maintaining the hESc phenotype and so it is important that the hESc are not overexposed to the Trypsin:EDTA. The cells should remain in clumps of approximately 5-10 cells to permit the formation of new hESc colonies.**

14. Transfer the hESc suspension to a sterile 15ml polystyrene centrifuge tube and centrifuge at 720 x g for 3 minutes at room temperature. Resuspend the hESc in 12ml of hESc medium.

15. Remove the MEF medium from the wells of the 6 well culture dish and add 2ml of the hESc cell suspension to each well. Transfer the culture dish to the cell culture incubator.
Tip: We frequently keep 1×6 well cell culture dish of hESc in continuous culture on an iMEF feeder layer. When confluent 1 x well is passaged to another 6 well culture dish with iMEFs, 2 x wells are taken to establish feeder-free cultures and 3 wells are cryopreserved in order to maintain low passage stocks of hESc.

Table 3: Formulation of hESc Feeder-Free Medium

hESc Feeder-Free Medium (500ml)	
DMEM:F12	480ml
BSA	0.5g
NEAA	5ml
B27 Supplement	10ml
N2 Supplement	5ml
β-Mercaptoethanol	91.7μl
Pass the medium through a 0.22μm filter, aliquot in 40ml volumes and store at -20°C	
Growth Factor Supplementation of Aliquots (40ml)	
100μg/ml (wt/vol) stock FGF2.	16μl
10μg/ml (wt/vol) stock Activin-A	40μl
10μg/ml (wt/vol) stock NT4	8μl
Store complete aliquots of hESc Medium at 4°C for up to 14 days	

Tip: Prior to assembling the components of the hESc Feeder-Free Medium transfer approximately 20ml of DMEM:F12 into a sterile 50ml centrifuge tube containing the 0.5g of BSA. Vortex the centrifuge tube well to promote complete dissolution of the BSA.

Feeder-Free hESc Culture:

1. Prepare 35mm and 6 well culture dishes by transferring 1ml of 50μg/ml fibronectin to each well and incubating for at least 16 hours at 4°C.
2. Remove and discard the hESc medium from 2 x wells of hESc iMEF cultures and rinse twice with DPBS.
3. Add 500μl of Trypsin:EDTA to each of the 2 wells and transfer to the cell culture incubator.
4. Check the hESc cultures every 2-3 minutes as if passaging the hESc cultures on to iMEF feeder cell layers. The Trypsin:EDTA can be left longer to permit more thorough dissociation of the hESc.
5. Transfer the cell suspension to a sterile 15ml polystyrene centrifuge tube and centrifuge at 720 x g for 3 minutes at room temperature.

6. Discard the supernatant and flick the bottom of the tube to disrupt the cell pellet. Resuspend the cells in a 2ml of hESc Feeder-Free Medium (Table 3).

7. Remove the 50µg/ml fibronectin from the 35mm culture dish and transfer to sterile 50ml centrifuge tube. **Tip: The 50µg/ml working solution of fibronectin can be reused up to three times if it remains at 4°C.**

8. Add the 2ml volume of hESc cell suspension to the well coated with fibronectin and check under the microscope. **Critical: To establish feeder-free cultures hESc should be kept at a high cell density transferring 2 x wells of a 6 well culture dish to 1×35mm culture dish to account for the inclusion of iMEFs. Cell-cell contact is important to maintain hESc phenotype, which is achieved through high cell density. Groups of 2-3 hESc not separated by the Trypsin:EDTA should still be present, however if large numbers of hESc remain aggregated together this will lead to the formation of embryoid bodies and differentiation.**

9. Return the 35mm culture dish to the cell culture incubator.
10. Check the hESc cultures every 24 hours under the microscope for adherence to the fibronectin-coated plastic and replace the hESc Feeder-Free Medium every 24-48 hours. **Tip: Feeder-free hESc adhere to the tissue culture plastic as colonies with prominent cell-cell contact. hESc appear larger and flatter than those cultured on iMEF feeder cell layers. Individual cells are seen as having characteristic large nucleus and prominent nucleoli. iMEF cells are often seen between hESc colonies.**

11. The feeder-free culture should be passaged at confluence after approximately 3-4 days. Remove the hESc Feeder-Free Medium from the well and add 500µl of pre-warmed Trypsin:EDTA. Incubate at room temperature and rock the 35mm culture dish back and forth gently. As soon as the hESc detach add 2ml of hESc Feeder-Free Medium to the well, draw up the hESc suspension and transfer to a sterile 15ml polystyrene centrifuge tube. Add another 2ml of hESc Feeder-Free Medium to the culture dish and rinse out the remaining cells from the well. Pool this with the hESc suspension in the 15ml tube. Centrifuge the hESc suspension at 720 x g for 3 minutes at room temperature and discard the supernatant. Flick the bottom of the centrifuge tube to disrupt the cell pellet and resuspend the hESc in 4ml volume of hESc Feeder-Free Medium. Transfer 2ml each to 2 x well of a 6 well cell culture dish coated with 50µg/ml fibronectin as described in step 1. Return the 6 well cell culture dish to the cell culture incubator. **Tip: It is important to maintain a high cell density as the hESc culture becomes established, plate cells at a density of approximately 0.8-1×10⁵cells/cm². Do not be tempted to passage the feeder-free hESc too early.**

12. Repeat steps 10-11 to expand the feeder-free hESc culture. The hESc become adapted to feeder-free culture with increasing passage and as such can be passaged at higher ratios. We routinely passage feeder-free hESc cultures at ratios of 1:3–1:4.

13. To cryopreserve feeder-free hESc cultures detach the cells as if passaging and resuspend the cell pellet in cryopreservation medium. Use 2ml of cryopreservation medium per 1 x confluent well of a 6-well dish (approximately 2×10⁶ cells). Aliquot 1ml volumes into pre-labelled cryovials, place the cryovials into the “Mr Frosty” and transfer to -80°C. After 24 hours transfer the cryovials to a liquid nitrogen cell storage dewar.

Table 4: Formulation of Directed Differentiation Basal Medium

Directed Differentiation Basal Medium (500ml)	
DMEM:F12	480ml
NEAA	5ml
B27 Supplement	10ml
ITS	5ml
β -Mercaptoethanol	917 μ l
Pass the medium through a 0.22 μ m filter, aliquot in 40ml volumes and store at -20°C. Thawed aliquots of Directed Differentiation Medium are stored at 4°C for up to 14 days	

Directed Differentiation Protocol

The directed differentiation protocol is carried out on feeder-free hESc which have been cultured off feeders for at least 3 passages to exclude contaminating iMEFs.

1. Coat 1×35mm culture dish by adding 1ml of 50 μ g/ml fibronectin and incubating overnight at 4°C.
2. Detach feeder-free hESc from the cell culture plastic using pre-warmed Trypsin:EDTA, centrifuge at 720 x g for 3 minutes at room temperature and resuspend to a cell density of 5×10⁵ cells/ml in hESc Feeder-Free Medium.
3. Remove the fibronectin solution from the well of the 35mm culture dish and add a 2ml volume of feeder-free hESc suspension.
4. Return the culture dish to the cell culture incubator.
5. After 24 hours check the hESc culture under the microscope. DDP should be initiated when the cultures are at approximately 80% confluence. Denote this as Day 1.
6. Prepare Day 1 medium by adding 6.25 μ l of 10 μ g/ml mouse recombinant Wnt3a and 12.5 μ l of 10 μ g/ml human recombinant Activin-A to 2.5ml of Directed Differentiation Basal Medium (Table 4).
7. Remove the hESc Feeder-Free Medium, rinse the feeder-free hESc culture three times with DPBS and add 2.5ml of Day 1 medium. Return to the cell culture incubator.
8. Day 2: add 6.25 μ l of 10 μ g/ml mouse recombinant Wnt3a; 6.25 μ l of 10 μ g/ml human recombinant Activin-A and 0.5 μ l of 100 μ g/ml human recombinant FGF2 to 2.5ml of Directed Differentiation Basal Medium.
9. Remove the Day 1 medium from the cell culture and replace with 2.5ml of Day 2 medium. Return to the cell culture incubator.
10. Day 3: add 6.25 μ l of 10 μ g/ml mouse recombinant Wnt3a; 2.5 μ l of 10 μ g/ml human recombinant Activin-A; 0.5 μ l of 100 μ g/ml human recombinant FGF2 and 10 μ l of 10 μ g/ml human recombinant BMP4 to 2.5ml of Directed Differentiation Basal Medium.
11. Remove the Day 2 medium from the cell culture and replace with 2.5ml of Day 3 medium. Return to the cell culture incubator.
12. Day 4: Analyse Stage 1 cultures if appropriate.

13. Add 0.5µl of 100µg/ml human recombinant FGF2; 10µl of 10µg/ml human recombinant BMP4; 25µl of 10µg/ml human recombinant follistatin and 0.5µl of 10µg/ml human recombinant NT4 to 2.5ml of Directed Differentiation Basal Medium.

14. Remove the Day 3 medium from the cell culture and replace with 2.5ml of Day 4 medium. Return to the cell culture incubator.

15. Transfer 1ml of 50µg/ml fibronectin to each of 5 x wells of a 6 well culture dish and place at 4°C.

16. Day 5: add 2µl of 100µg/ml human recombinant FGF2; 40µl of 10µg/ml human recombinant BMP4; 100µl of 10µg/ml human recombinant follistatin and 2µl of 10µg/ml human recombinant NT4 to 10ml of Directed Differentiation Basal Medium.

17. Remove the Day 4 medium from the cell culture and add 500µl of pre-warmed Trypsin:EDTA. Incubate at room temperature and rock the 35mm culture dish back and forth gently. As soon as the cells detach add 2ml of Directed Differentiation Basal Medium to the well, draw up the cell suspension and transfer to a sterile 15ml polystyrene centrifuge tube. Add another 2ml volume of Directed Differentiation Basal Medium to the culture dish and rinse out the remaining cells from the well. Pool this with the cell suspension. Centrifuge the cell suspension at 720 x g for 3 minutes at room temperature and discard the supernatant. Flick the bottom of the centrifuge tube to disrupt the cell pellet and resuspend the cells in 10ml of Day 5 Medium. Remove the 50µg/ml fibronectin solution and transfer 2ml of the cell suspension to each of the 5 wells. Return the culture dishes to the cell culture incubator.

18. Day 6: add 2µl of 100µg/ml human recombinant FGF2; 40µl of 10µg/ml human recombinant BMP4; 100µl of 10µg/ml human recombinant follistatin and 2µl of 10µg/ml human recombinant NT4 to 10ml of Directed Differentiation Basal Medium.

19. Remove the Day 5 medium from the 5 cell culture dishes and replace each with 2ml of Day 6 medium. Return to the cell culture incubator.

20. Day 7: add 2µl of 100µg/ml human recombinant FGF2; 40µl of 10µg/ml human recombinant BMP4; 100µl of 10µg/ml human recombinant follistatin and 2µl of 10µg/ml human recombinant NT4 to 10ml of Directed Differentiation Basal Medium.

21. Remove the Day 6 medium from the 5 cell culture dishes and replace each with 2ml of Day 7 medium. Return to the cell culture incubator.

22. Transfer 10ml of 50µg/ml fibronectin and 10ml of 0.1% (vol/vol) gelatin to a sterile 50ml polystyrene centrifuge tube and mix well by pipetting up and down.

23. Transfer 1ml of the mixed fibronectin:gelatin solution to 20 x wells of 6 well culture dishes and place at 4°C.

24. Day 8: add 8µl of 100µg/ml human recombinant FGF2; 160µl of 10µg/ml human recombinant BMP4 and 8µl of 10µg/ml human recombinant NT4 to 40ml of Directed Differentiation Basal Medium.

25. Remove the Day 7 medium from the cell cultures and add 500µl of pre-warmed Trypsin:EDTA to each of the 5 wells. Incubate at room temperature; rocking the cell culture dish back and forth gently. As soon as the cells detach add 2ml of Directed Differentiation Basal Medium to each cell culture dish, draw up the cell suspension and transfer to a sterile 15ml polystyrene centrifuge tube. Add another 2ml

of Directed Differentiation Basal Medium to the culture dishes and rinse out the remaining cells from the well. Pool this with the cell suspension. Centrifuge the cell suspension at 720 x g for 3 minutes at room temperature and discard the supernatant. Flick the bottom of the centrifuge tube to disrupt the cell pellet and resuspend the cells in 40ml of Day 8 Medium. Remove the fibronectin:gelatin solution and transfer 2ml of the cell suspension to each of the 20 wells of the culture dishes. Return the culture dishes to the cell culture incubator.

26. Day 9: Analyse Stage 2 cultures if appropriate.

27. Add 8µl of 100µg/ml human recombinant FGF2; 80µl of 10µg/ml human recombinant BMP4; 80µl of 10µg/ml human recombinant GDF5 and 8µl of 10µg/ml human recombinant NT4 to 40ml of Directed Differentiation Basal Medium.

28. Remove the Day 8 medium from the cell cultures and replace with 2ml of Day 9 medium to each well. Return to the cell culture incubator.

29. Day 10: add 8µl of 100µg/ml human recombinant FGF2; 80µl of 10µg/ml human recombinant BMP4; 80µl of 10µg/ml human recombinant GDF5 and 8µl of 10µg/ml human recombinant NT4 to 40ml of Directed Differentiation Basal Medium.

30. Remove the Day 9 medium from the cell cultures and replace with 2ml of Day 10 medium to each of the wells. Return to the cell culture incubator.

31. Day 11: add 8µl of 100µg/ml human recombinant FGF2; 160µl of 10µg/ml human recombinant GDF5 and 8µl of 10µg/ml human recombinant NT4 to 40ml of Directed Differentiation Basal Medium.

32. Remove the Day 10 medium from the cell cultures and replace with 2ml of Day 11 medium to each well. Return to the cell culture incubator.

33. Transfer 1ml of 0.1% (vol/vol) gelatin solution to 30 x wells of 6 well cell culture dishes and place at 4°C.

34. Day 12: add 12µl of 100µg/ml human recombinant FGF2; 240µl of 10µg/ml human recombinant GDF5 and 12µl of 10µg/ml human recombinant NT4 to 30ml of Directed Differentiation Basal Medium.

Note: This makes a 2 X concentration of Day 12 medium.

35. Remove the Day 11 medium from the cell cultures and add 500µl of pre-warmed Trypsin:EDTA to each well. Incubate at room temperature; rocking the cell culture dishes back and forth gently. As soon as the cells detach add a 2ml volume of Directed Differentiation Basal Medium to each cell culture dish, draw up the cell suspension and transfer in equal volumes to 2 x sterile 50ml polystyrene centrifuge tubes. Add another 2ml volume of Directed Differentiation Basal Medium to the culture dishes and rinse out the remaining cells from the well. Transfer this cell suspension in equal volumes to the 2x50ml centrifuge tubes. Centrifuge the cell suspension at 720 x g for 3 minutes at room temperature and discard the supernatant. Flick the bottom of the centrifuge tube to disrupt the cell pellet and resuspend the cells in 30ml of Directed Differentiation Basal Medium. Remove the 0.1%(vol/vol) gelatin solution and transfer 1ml of the cell suspension to each of the 30 wells of the 6 well culture dishes.

36. Add 1ml of the 2 X Day 12 medium to each of the 30 wells. Return the culture dishes to the cell culture incubator.

37. Day 13: add 12µl of 100µg/ml human recombinant FGF2; 240µl of 10µg/ml human recombinant

GDF5 and 12µl of 10µg/ml human recombinant NT4 to 30ml of Directed Differentiation Basal Medium.
Note: This makes a 2 X concentration of Day 13 medium.

38. Remove the Day 12 medium from the cell cultures and replace with 1ml each of Directed Differentiation Basal Medium to each of the 30 wells of the 6 well cell culture dishes

39. Add 1ml each of the 2 X Day 13 medium to each of the 30 wells. Return to the cell culture incubator.

40. Day 14 terminate the cultures and carry out Stage 3 (end point) analyses.

REFERENCES

A chemically-defined protocol for generating chondrocytes from human embryonic stem cells
[Rachel Oldershaw](#) , [Melissa Baxter](#) , [Emma Lowe](#) , [Nicola Bates](#) , [Lisa Grady](#) , [Francesca Soncin](#) ,
[Daniel Brison](#) [Timothy Hardingham](#) & [Susan Kimber](#)
Protocol Exchange (2010) Published online 30 November 2010